


# FLoid™ Cell Imaging Station quick start guide

## What's in the box?


FLoid™ cell imaging station  
(Cat. No. 4471136)  
Power cord  
Keyboard  
Mouse  
USB flash drive  
(w/ User manual)  
Quick start guide  
Printer (optional)



## 1. Connect accessories

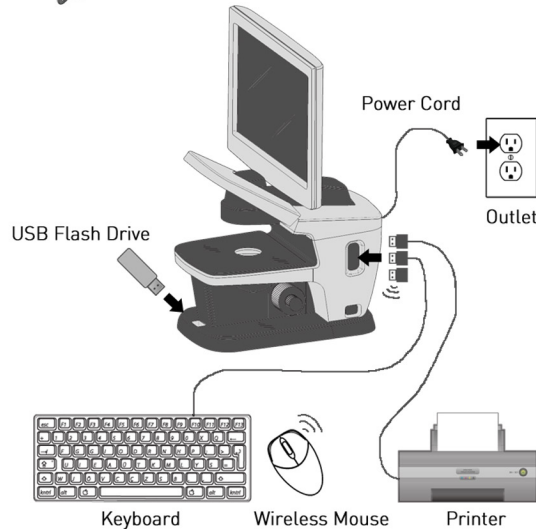

## 2. Power on


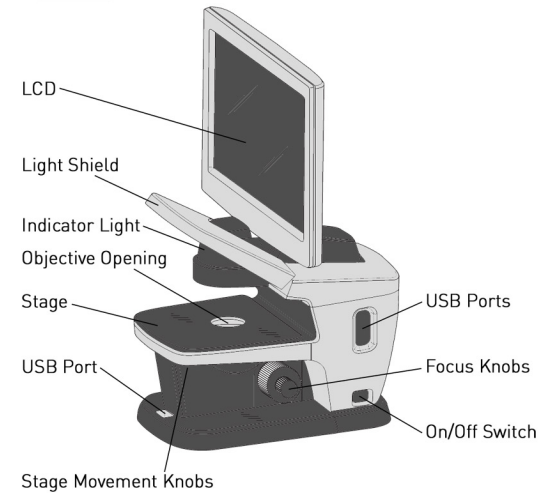


# FLoid™ Cell Imaging Station quick start guide



## 3. Capture Image

- Place sample on stage over objective opening
- Turn on light by selecting light bulb icon
- Focus image using focus knobs
- Capture image(s) then press "Processing" button

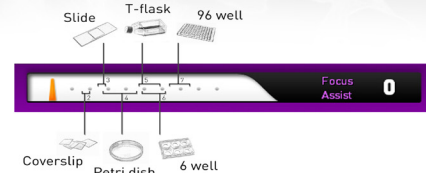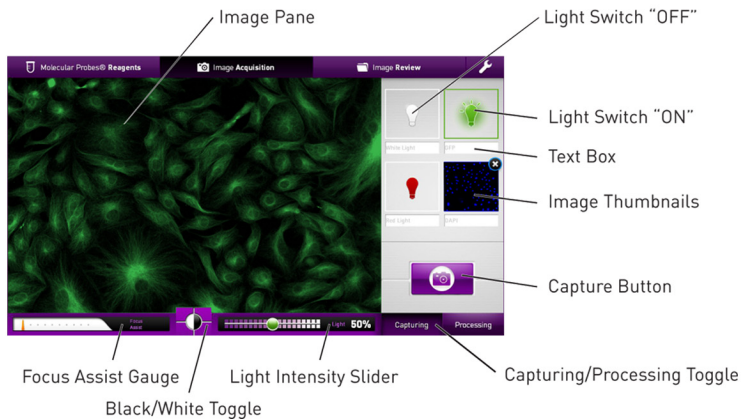

## 4. Process Image

- Adjust brightness and contrast as needed
- Save image to USB flash drive and print image

